



	<u>Sample</u>	<u>40 U/ml tPA</u>	<u>PAI-1 Assay Buffer</u>
Sample	20 µl	20 µl	--
0 U/ml PAI-1	--	20 µl	20 µl
40 U/ml PAI-1	--	--	40 µl

Notes:

- Citrate treated platelet poor plasma must be used for the assay. Platelet contamination may cause spurious results.
 - PAI-1 is unstable and must be processed and frozen within 3 hr of specimen collection.
 - There is large diurnal variation of PAI-1 activity in plasma, which should be taken into consideration when designing clinical studies and routine applications. It is recommended that specimens should be the early morning fasting specimen.
 - If PAI-1 activity is above 40 U/ml, dilute the sample(s) with PAI-1 Assay Buffer and mark the dilution factor.
- 3. Acidification & Neutralization Step:** Add 40 µl of PAI-1 Acidify Buffer to all three eppendorf tubes including **Sample, 0 U/ml PAI-1** and **40 U/ml PAI-1** (from Step 2). Mix well and incubate at 37°C for 20 min, protected from light. Add 80 µl of PAI-1 Assay Buffer to all the three tubes and mix well. The "**Sample**" tube is now ready for the assay. Add 10 µl of the Sample into two parallel wells of a clear 96-well half area plate designed as "**PAI-1 Sample**" and "**PAI-1 Sample Background Control**".

Note: Equilibrate the clear 96-well half area plate to 37°C before adding the Sample(s).

4. Standard Curve Preparation:

- Intermediate "PAI-1" Standard Preparation:** Prepare various intermediate PAI-1 Standards including 10 U/ml, 20 U/ml and 30 U/ml using PAI-1 Acidify Buffer treated 0 U/ml PAI-1 and 40 U/ml PAI-1 (from Step 3) according to the table below. Mix well.

	<u>0 U/ml PAI-1</u>	<u>40 U/ml PAI-1</u>
10 U/ml PAI-1	45 µl	15 µl
20 U/ml PAI-1	30 µl	30 µl
30 U/ml PAI-1	15 µl	45 µl

- Add 10 µl of 0 U/ml PAI-1, 10 U/ml PAI-1, 20 U/ml PAI-1, 30 U/ml PAI-1 and 40 U/ml PAI-1 Standards (from Step 3 & 4) into wells of clear 96-well half area plate.

- 5. Reaction Mix Preparation:** Prepare a 25-fold dilution of the Plasminogen stock solution with PAI-1 Assay Buffer (i.e. add 4 µl of Plasminogen stock solution with 96 µl PAI-1 Assay Buffer). Prepare 10-fold dilution of the reconstituted Substrate Mix with PAI-1 Assay Buffer (i.e. add 10 µl of reconstituted Substrate Mix with 90 µl PAI-1 Assay Buffer).

Prepare Reaction Mix (for both Standard(s) & PAI-1 Sample wells) and Background Mix (for PAI-1 Sample Background Control wells) according to the table below. Make sufficient amount of each type of mix to add 40 µl to all assay wells of that type.

	<u>Reaction Mix</u>	<u>Background Mix</u>
Diluted Plasminogen	10 µl	-- µl
Diluted Substrate Mix	10 µl	10 µl
PAI-1 Assay Buffer	20 µl	30 µl

Mix well. Add 40 µl of Reaction Mix to PAI-1 Standard(s) & PAI-1 Sample wells and 40 µl of Background Mix to PAI-1 Sample Background Control well(s). Mix well and incubate at 37°C for 90 min, protected from light. The total volume of each well is 50 µl. After 90-min incubation, add 50 µl of PAI-1 Stop Buffer to all wells containing PAI-1 Sample(s), PAI-1 Sample Background Control and PAI-1 Standards. Mix well.

Note: Equilibrate the PAI-1 Stop Buffer to 37°C before adding to the wells.


- 6. Measurement:** Measure the colorimetric signal (OD 405 nm) at 37°C in end-point mode.

- 7. Calculation:** Subtract the "40 U/ml PAI-1" Standard reading from all Standards readings. Plot the PAI-1 Standard Curve. Subtract the PAI-1 Sample Background Control reading from all PAI-1 Sample readings to get the corrected PAI-1 Sample readings. Apply the corrected PAI-1 Sample readings to the PAI-1 Standard Curve to obtain the corresponding PAI-1 activity (U/ml) as:

$$\text{Sample PAI-1 Activity} = B * D = \text{U/ml}$$

Where: **B** = PAI-1 activity from the Standard Curve (U/ml)
D = Sample dilution factor (D=1 for undiluted Sample(s))

Unit Definition: One unit of PAI-1 activity was defined as the amount of PAI-1 that inhibits one unit of tPA activity under Assay conditions.

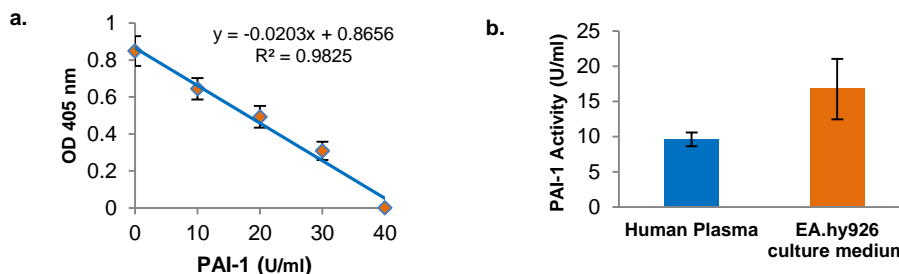

Figures: a. PAI-1 Standard Curve. b. Measurement of PAI-1 Activity in pooled human plasma (citrate treated platelet poor plasma) and EA.hy926 culture medium. All assays were performed following kit protocols.

VIII. Related Products:

Tissue Plasminogen Activator Activity Assay Kit (K178)
Urokinase Activity Fluorometric Assay Kit (K728)

Plasmin Activity Assay Kit (Fluorometric) (K381)
Plasmin Activity Assay Kit (Colorimetric) (K945)

FOR RESEARCH USE ONLY! Not to be used on humans