



EZClick™ Myristoylated Protein Assay Kit (FACS/Microscopy)

rev 07/20

(Catalog # K497-100; 100 assays; Store at -20°C)

I. Introduction:

Myristoylation is a co- and post-translational modification of proteins. Myristoylated proteins are involved in cellular trafficking, cellular functions and oncogenesis. In this modification, Myristic acid, a saturated fatty acid is covalently bond to a protein through the action of *N*-myristoyl transferase (NMT). NMT transfers the myristate moiety from Myristoyl-CoA to an *N*-terminal glycine of a target protein. NMT is associated with carcinogenesis. The upregulation of its expression and activity is observed in colorectal cancer, gall-bladder cancer and brain tumors. BioVision offers **EZClickTM Myristoylated Protein Assay Kit**, a highly specific, simple and robust method for labeling and detection of myristoylated proteins. We use Alkynyl Myristic Acid that is fed directly into the cells and gets incorporated into proteins during or post translation, followed by click reaction with an azide-containing dye. This system offers a powerful method for imaging localization, trafficking, and dynamics of myristoylated proteins, or detection by FACS for quantitative studies. Labeled myristoylated proteins can be appropriate excitation sources, or enriched by immunoprecipitation with biotin-azide or antibodies prior to proteomic analysis. We provide sufficient materials for 100 assays in a 96-well plate format.

II. Applications:

- Identification and localization of myristic acid modified proteins.
- · Detection and quantification of biosynthesis, dynamics and turnover of myristic acid modified proteins
- Screening for genotoxic compounds and effectors of protein modifications

III. Sample Type:

• Suspension or adherent cells

IV. Kit Contents:

Components	K497-100	Cap Code	Part Number
EZClick [™] Wash Buffer (10X) Fixative Solution	25 ml 10 ml	NM WM	K497-100-1 K497-100-2
Permeabilization Buffer (10X) EZClick [™] Myristic Acid Label (1000X)	25 ml 10 μl	NM/Blue White	K497-100-3 K497-100-4
Copper Reagent (100X)	100 µl	Blue	K497-100-5
EZClick [™] Fluorescent Azide (100X)	100 µl	Green/Amber	K497-100-6
Reducing Agent (20X)	500 µl	Yellow	K497-100-7
EZClick [™] Total DNA Stain (1000X)	10 µl	Blue/Amber	K497-100-8

V. User Supplied Reagents and Equipment:

- Tissue culture vessels and appropriate culturing media
- Phosphate Buffered Saline (PBS, pH 7.4)
- Sterile 0.1% Gelatin Solution (optional, only required for suspension cells)
- Flow cytometer equipped with laser capable of excitation at 488 nm wavelength (FL-1)
- · Fluorescence microscope capable of excitation and emission at 440/490 nm and UV filter

VI. Storage Conditions and Reagent Preparation:

Upon arrival, store the entire kit at -20°C protected from light. Briefly centrifuge small vials prior to opening. Read the entire protocol before performing the assay.

- 10X Wash Buffer and 10X Permeabilization Buffers: Thaw at 37°C to dissolve completely. Dilute the 10X stocks 1:10 in sterile water, mix well. Store at 4°C.
- Fixative Solution: Divide into aliquots and store at -20°C, protected from light.
- Remaining components: Store at -20°C protected from light. While in use, keep on ice and minimize light exposure.

VII. Myristoylated Protein Assay Protocol:

Notes:

This assay was developed with HeLa (adherent) and Jurkat (suspension) cells and can be modified for any cell line. The protocol below refers to a 96-well tissue culture plate. Adjust volumes accordingly for other plate formats. The assay volume is 100 µl. Growth conditions, cell number per well and other factors may affect the incorporation rate of the Protein Label. Therefore optimize the assay for your cell type. We suggest an initial test of several EZClick[™] Myristic Acid Label concentrations to find best conditions for your experimental design. Avoid stressing the cells by washes or temperature changes prior to incubation with EZClick[™] Myristic Acid Label. All steps should be carried out at room temperature (RT) unless otherwise specified; equilibrate all buffers to RT prior to the experiment

1. Labeling of control and experimental cells: method with drug pre-incubation:

- a. Obtain cell suspension of desired density and seed directly into tissue culture vessels or on coverslips for high resolution microscopy. <u>To immobilize suspension cells for microscopy</u>: add 100 µl of 0.1% gelatin solution directly into the wells. Tilt the plate to cover the entire well surface and place it in a tissue culture hood for 1 hour. Gently remove the gelatin solution and seed your cells. Allow the cells to recover overnight before the treatment. Next day, treat the cells with appropriate effectors according to your protocol; *do not add treatment to the positive and negative control cells*. <u>Negative Control Cells</u> (Unstained Cells): cells not exposed to Myristic Acid Label or EZClick[™] Fluorescent Azide), <u>Background Control Cells</u> (Cells are not exposed to the EZClick[™] Myristic Acid Label, EZClick[™] Reaction only), <u>Positive Control Cells</u> (Cells are incubated with 1X EZClick[™] Myristic Acid Label and EZClick[™] Reaction).
- b. Next day, for adherent cells, remove the media directly. For suspension cells, centrifuge cells at 500 x g for 5 min, and discard the supernatant. Then replace it with fresh aliquots containing EZClick[™] Myristic Acid Label (1000X) diluted to 1X final concentration with





culture medium and add into the experimental and positive control cells respectively. *Do not add* the EZClick[™] Myristic Acid Label into the **Negative Control Cells**.

- c. Add treatments and incubate the cells for additional 1 day or for the period of time required by your experimental protocol. Do not remove the drug containing media while incubating with 1X EZClick[™] Myristic Acid Label to avoid potential reversibility of drug action on label incorporation.
- d. To terminate the experiment, For adherent cells: Remove the media and rinse the cells once with 100 µl of 1X PBS, discard the supernatant. For suspension cells: Centrifuge the cells at 500 x g for 5 min to deposit the cells onto the surface. Tilt the plate and gently remove the media with a pipette tip. It is important to avoid excessive centrifugation speeds, which can damage the cells. Make note of the place that is used, and perform subsequent aspirations from the same place. Pellet the suspension cells at 500 x g for 5 min throughout the entire protocol!

2. Fixation and Permeabilization:

- a. For adherent cells: Add 100 µl of Fixative Solution to each well and incubate the cells for 15 min at RT protected from light. Remove the fixative and wash the cells once with 100 µl of 1X Wash Buffer, remove the wash. Add 100 µl of 1X Permeabilization Buffer and incubate the cells for 10 min at RT. Remove the supernatant. Proceed to EZClick[™] Myristic Acid reaction and total DNA staining.
- b. For suspension cells: Re-suspend the cells in 100 µl of Fixative Solution and incubate for 15 min at RT protected from light. Centrifuge the cells at 500 x g for 5 min, remove the fixative and wash the cells once with 100 µl of 1X Wash Buffer. Centrifuge the cells at 500 x g for 5 min, discard the supernatant and re-suspend the cells in 100 µl of 1X Permeabilization Buffer. Incubate the cells for 10 min at RT. Centrifuge the cells at 500 x g for 5 minutes, remove the supernatant. Proceed to EZClick™ Myristic Acid reaction and total DNA staining.

3. EZClick™ Myristic Acid reaction and total DNA staining:

a. Reaction Cocktail: Prepare 1X EZClick[™] Reaction Cocktail according to the table below. Volumes should be multiplied by number of samples and reagents added in the exact order. Use the Reaction Cocktail within 15 min of preparation. *Cells should be protected from light during, and following the EZClick[™] reaction and DNA staining.*

	Amount per Reaction
PBS	93 µl
Copper Reagent (100X)	1 µl
EZClick [™] Fluorescent Azide (100X)	1 µl
Reducing Agent (20X)	5 µl

- b. EZClick™ Myristic Acid Reaction: For Negative Control Cells, add 100 µl of 1X PBS. For Background Control Cells, Positive Control Cells and Experimental Cells: Add 100 µl of 1X EZClick™ Reaction Cocktail to each sample and incubate the cells for 30 min at RT protected from light. Remove the Reaction Cocktail and wash cells three times in 100 µl of 1X Wash Buffer. Remove the 1X wash and suspend the cells in 100 µl of 1X PBS. Proceed to DNA staining. If no DNA staining is desired, proceed to Microscopic or FACS analysis. DNA staining: Prepare 1X dilution of Total DNA Stain and add 100 µl per well. Incubate the cells for 20 min at RT, or refrigerate at 4°C protected from light. Remove the DNA stain solution; wash the cells once with 100 µl PBS. Note: cells are compatible with all methods of slide preparation including wet mount or prepared mounting media.
- 4. Fluorescence Microscope Imaging: Analyze samples for green fluorescence generated by EZClick[™] labeled Myristic Acid and for blue fluorescence by nuclear DNA. FACS analysis: Harvest cells by preferred method and wash with 0.5 ml of ice-cold PBS. Resuspend pellets with 100 µl of ice-cold PBS and analyze samples for green fluorescence generated by EZClick[™] Myristic Acid addition during EZClick[™] reaction.



CL 05025 110 + 170 (400) 402 1000 TE (400) 402 1001 1

Figures Analysis of metabolic labeling of Myristic Acid on proliferating cells. HeLa (10⁵ cells/ ml) and Jurkat (1X10⁶ cells/ml) cells respectively were incubated overnight with fresh aliquots of media containing EZClickTM Myristic Acid Label. Cells were then processed and analyzed by Microscopy and FACS according to the kit protocol. (A) HeLa cells: Upper panel corresponds to the Azide only Background fluorescence. The lower panel shows that the green fluorescence corresponds to the EZclickTM Myristic Acid labeling. Nuclear staining in both panels confirms that green signal is a result of Myristic Acid Label incorporation. (B) Jurkat cells: FACS analysis of negative control (Black), background (EZClick only, Green), positive control (Myristic Acid Label and EZClick, Pink). Signal measured in FL-1 channel clearly shows the Myristic Acid Labeling of Protein.

VIII. Related EZClick[™] Products:

Global RNA Assay Kit (K718) EdU Cell Proliferation/DNA Synthesis Kit (K946) Global Phospholipid Assay Kit (K717)

.. ..

.

FOR RESEARCH USE ONLY! Not to be used on humans.